



Kleenpak™ Capsules with Ultipor® N66 Membrane

Description

Pall Kleenpak sterilizing-grade capsule filters with intrinsically hydrophilic Ultipor N66 membrane are self-contained sanitary filters designed for convenient small-batch sterile filtration of most **biopharmaceutical** solutions. Rugged, compact design ensures reliability.



Features and Benefits

- ▶ 100% integrity and pressure tested
- ▶ 0.2 micron NFP sterilizing-grade
- ▶ 0.1 micron NTP sterilizing-grade for bacteria and mycoplasma removal
- ▶ Low filter extractables
- ▶ Excellent solvent compatibility
- ▶ Sanitary vent and drain valves with non-twist hose barbs
- ▶ Low hold-up volumes (< 1 mL)
- ▶ Gamma-irradiatable, autoclavable, and pre-sterilized versions available for maximum flexibility
- ▶ Autoclavable versions can be repeatedly autoclaved
- ▶ Manufactured for use in conformance with cGMP
- ▶ ISO 9000 certified Quality system
- ▶ Pharmaceutical P optimized
- ▶ Certificate of test provided

Quality and Bio-Safety Biological Tests

Integrity

- ▶ Every filter tested during manufacture. Test correlated to microbial retention

Biological Tests

- ▶ Meets USP Biological Reactivity test, in vivo, for Class VI-121 °C Plastics

Effluent Quality Tests*

- ▶ Meets Cleanliness per USP Particulates in Injectables
- ▶ Non-Fiber-Releasing
- ▶ Non-Pyrogenic per USP Bacterial Endotoxins (< 0.25 EU/mL)
- ▶ Meets Total Organic Carbon and Water Conductivity per USP Purified Water; pH per Sterile Purified Water

Autoclave Resistance

- ▶ Lot samples multi-cycle autoclave challenged

* Per lot sample or rinse-flush aliquots.

Specifications

Materials of Construction

Membrane	Nylon 6,6
Support and Drainage	Polyester
Core and Cage	Autoclavable Version: Polypropylene G and S Grades: Polyester
End Caps and Shell	Polyester
Vent and Drain Valve O-ring	Ethylene Propylene (EPDM)

Microbial Removal Ratings¹

NF 0.2 µm, sterilizing-grade, *Brevundimonas diminuta*

NT 0.1 µm, sterilizing-grade,²*Brevundimonas diminuta*

¹ Assembly lot samples retain > 10⁷ cfu/cm² of *Brevundimonas diminuta* per modified ASTM F838-83 and FDA guidelines.

² Also retains > 10⁷ cfu/cm² *Acholeplasma laidlawii* mycoplasma.

Nominal Filter Areas

Size Code	KA1	KA2	KA3	KA4
NF	0.05 m ² (0.53 ft ²)	0.1 m ² (1 ft ²)	0.2 m ² (2.1 ft ²)	0.5 m ² (5.3 ft ²)
NT	0.045 m ² (0.4 ft ²)	0.08 m ² (0.8 ft ²)	0.2 m ² (2.1 ft ²)	0.4 m ² (4 ft ²)

Nominal Dimensions

Size Code	KA1	KA2	KA3	KA4
Maximum Diameter of Bowl (including Valves)	94 mm (3.7 in.)	94 mm (3.7 in.)	109 mm (4.2 in.)	109 mm (4.2 in.)
Length (including Sanitary Connection* [Code 1 and 8])	117 mm (4.6 in.)	158 mm (6.2 in.)	174 mm (6.8 in.)	286 mm (11.2 in.)


* Sanitary connection is Tri-Clamp compatible. Tri-Clamp is a trademark of Alfa Laval, Inc

Operating Conditions³

Maximum Pressure and Temperature	5.2 barg (75 psig) to 40 °C
Maximum Differential Pressure	4.1 barg (60 psid) at 40 °C

³ Using compatible liquids. Recommended maximum 3.5 barg (50 psig) in continuous air and gas service. KA*NT filters may be forward flow tested at 5171 mbar (75 psig) at ambient temperature.

Typical Liquid Flow Rates⁴

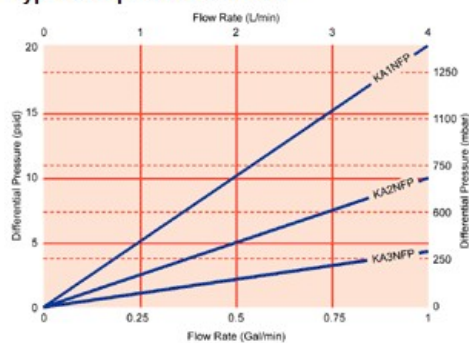

⁴ Typical initial clean ΔP; water at 20 °C, 1 cP. Values shown are for 38 mm (1½ in.) sanitary flange connections. Values with other connections are available on request. For assistance, contact your local Pall representative.

Aqueous Extractables (NVR)⁵

Size Code	KA1	KA2	KA3	KA4
Non irradiated	7 mg	10 mg	12 mg	20 mg
Pre-sterilized	6 mg	12 mg	22 mg	N/A

⁵ Water at 20 – 25 °C after autoclaving.
N/A = Filter not available

Steam Autoclaving⁶ (Non G and S Grades)

Cumulative Autoclave Time	16 hours at 125 °C 4 hours at 140 °C
---------------------------	---

⁶ Laboratory tests (1-hour cycles) establish multi-cycle. Filters should be qualified in actual use. Contact Pall for recommended procedures

Warning: Kleenpak filters must not be steam-sterilized in situ by passing steam through under pressure.

Sterilization

Gamma-irradiation Maximum Dosage (G Grade)	50 kGy
--	--------

Note: Contact Pall for specific part number availability.

S and G grades available in KA1 and KA2 with inlet and outlet options 1 and 2; KA3 with with inlet and outlet options 1 and 6.

Autoclavable grades available in KA1 and KA2 with inlet and outlet options 1 and 2; KA3 and KA4 with with inlet and outlet options 1 and 6.

Ordering Information

Ordering Information							
KA			P				
Code	Size and Area	Code	Removal Ratings	Code	Inlet and Outlet Connections	Code	Options
1	See above dimensions table	NF	0.2 µm sterilizing-grade	1	1 - 1½ in. sanitary flange	Blank	Autoclave service
2		NT	0.1 µm sterilizing-grade	2	6 - 13 mm (¼ - ½ in.) stepped hose barb	G	Unsterilized, Gamma-irradiatable
3				6	14 mm (½ in.) hose barb	S	Pre-sterilized
4				8	¾ - 1 in. fractional sanitary flange		

Option P8 available in KA2, autoclave service.

Contact Information

Pall Office(s)

Biopharm East
770 Pennsylvania Drive
Suite 100
Exton, PA 19341
USA
Phone: 610 458 9500
Fax: 610 458 9585

Biopharm West
1630 Industrial Park Street
Covina, CA 91722
Phone: 888.426.7255
Fax: 610.458.8090

World Headquarters
25 Harbor Park Drive
Port Washington, NY 11050
USA
Phone: (516) 484-3600
Alternate Phone: **1-800-289-7255**
Fax: (516) 801-9754
[Driving Directions](#)
[Map](#)

Biopharm Southeast
25 Harbor Park Drive
Port Washington, NY 11050
USA
Phone: 516.484.3600
Alternate Phone: 1.800.289.7255

Pall Canada Ltd. -Quebec
2535 De Miniac
Ville St. Laurent
Quebec, H4S 1E5
Phone: +1 514 332-7255
Alternate Phone: 800-435-6268 (Canada Only)
Fax: +1 514 332-0996
Alternate Fax: 800-808-6268 (Canada Only)

Distributor(s)

Enpro, Inc.
121 S. Lombard Road
Addison, IL 60101
Phone: 630-629-3504
Fax: 630-629-3512
Email: inquiry@enproinc.com
Website: <http://www.enproinc.com>

Enpro, Inc.
8888 Keystone Crossing
Suite 1300
Indianapolis, IN 46240
Phone: (317) 872-1422
Fax: (317) 872-1554
Email: inquiry@enproinc.com
Website: <http://www.enproinc.com>

Enpro, Inc.
7300 West 110th St., 7th Floor
Overland Park, KS 66210
USA
Phone: (913) 327-8400
Fax: (913) 327-8401
Email: inquiry@enproinc.com
Website: <http://www.enproinc.com>

Enpro, Inc.
7825 Washington Ave., Ste.
504

Enpro, Inc.
1001 Craig Road
Suite 260

Enpro, Inc.
11329 P St., Ste 110
Omaha, NE 68137

Bloomington, MN 55439
USA
Phone: (952) 943-3930
Fax: (952) 943-8022
Email: inquiry@enproinc.com
Website:
<http://www.enproinc.com>

St. Louis, MO 63146
USA
Phone: (314) 878-6161
Fax: (314) 878-0470
Email: inquiry@enproinc.com
Website:
<http://www.enproinc.com>

USA
Phone: (402) 597-2306
Fax: (402) 597-2320
Email: inquiry@enproinc.com
Website:
<http://www.enproinc.com>

Enpro, Inc.
3030 Northwest Expressway
Ste. 200, #121
Oklahoma City, OK 73112
USA
Phone: (405) 728-9500
Fax: (405) 728-9520
Email: info@enproinc.com
Website:
<http://www.enproinc.com>

Enpro, Inc.
5865 Ridgeway Center Pkwy.,
Ste. 300
Memphis, TN 38120-4004
USA
Phone: (901) 685-2929
Fax: (901) 685-2930
Email: inquiry@enproinc.com
Website:
<http://www.enproinc.com>

Enpro, Inc.
15950 N. Dallas Pkwy., Ste.
400
Dallas, TX 75248
USA
Phone: (972) 361-8320
Fax: (972) 361-8321
Email: inquiry@enproinc.com
Website:
<http://www.enproinc.com>

Inseco, Incorporated
Julia Industrial Park
A Street – Matadero Corner
San Juan, PUERTO RICO
00920
Tel: 787.781.2655
Fax: 787.793.3620
E-Mail:
Nestor_Guedez@insecopr.com
Contact: Nestor Guedez

© 2012, Pall Corporation, Pall, and other names are trademarks of Pall Corporation.
® indicates a registered trademark in the USA.